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Equipment Safety Instruction: Chemical storage cabinet

Task	Hazard	Risk	Control
Before storing chemicals	Chemical containers reuse	<ul style="list-style-type: none">- Generation of heat- Fire- Explosion- Contaminate	<ul style="list-style-type: none">- If reuse an empty bottle, deface the original label and create a new label- Keep a fire extinguisher handy- Wear the proper PPE
Chemical storage	Chemicals are stored in alphabetical order	<ul style="list-style-type: none">- Generation of heat- Fire- Explosion- Contaminate	<ul style="list-style-type: none">- Review the SDS- Separate and store chemicals by chemical hazard and compatibility group- All chemicals should be GHS labeled- Provide a secondary container
	Chemical spills	<ul style="list-style-type: none">- Skin burns or irritation caused by contact with the hazardous chemicals- Poisoning by absorption through the skin of a toxic substance	<ul style="list-style-type: none">- Wear the proper PPE- Avoid storing all chemicals above eye level and shoulder height- Preparation of a first aid kit and spill kit
	Chemical container's cap not tightly closed	<ul style="list-style-type: none">- Accidental inhalation or ingestion, vapors can also irritate the eyes	<ul style="list-style-type: none">- Wear the proper PPE- Keep all containers of chemicals closed when not in use
	Use of expired chemicals	<ul style="list-style-type: none">- Cancer occurring years after exposure to a carcinogenic substance	<ul style="list-style-type: none">- Use first-in, first-out system- Date all chemicals when received and when opened- Note expiration dates on chemicals

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