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Equipment Safety Instruction: Water bath

Task	Hazard	Risk	Control
Switch on equipment	Electricity	- Workers electrocute	- Use a properly grounded electrical outlet with the proper voltage and current handling capacity
		- Liquid spill cause an electric shock	- Place water bath on flat and level surface - Wear the proper PPE
		- Fire and explosion	- Do not store volatile flammable materials near a water bath
During the operation	Electricity	- Fire and explosion	- Do not leave a water bath unattended when in use
	Container is not completely close	- Liquid spillage - Damage to water bath	- Make sure water bath is filled to an appropriate level, Wear the proper PPE - Preparation of a spill kit
	Work with hazardous substances	- Injury through exposure to aerosols with chemical and biological hazards	- Wear the proper PPE - Request labels for water bath when in use
	Chipped or cracked glassware	- Injury through exposure to aerosols with chemical and biological hazards	- Inspect all glassware to be used in the centrifuge (chipped/cracked) - Preparation of a spill kit, first aid kit
	Thermal	- Worker injured	- If water temperature is above 50°C be very careful as skin contact with the water - If changing the temperature of the water bath, label the bath with the temperature you will be using it at and the date it will be used
	Add a tap water to the water bath	- Water bath is not operate at the desired temperature	- Filled with distilled water only

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Task	Hazard	Risk	Control
Maintenance		- Contaminate	- Preparing a cleaning schedule - Water baths should have the water changed in them regularly and wiped out with Virkon solution - Decontaminate by raising the temperature to 90°C or higher for 30 minutes once a week - Unplugging when not in use

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